

EnVision Multilabel Plate Reader



Is Your EnVision System Ready for AlphaPlex?

Alpha Technology is our versatile, homogeneous, easy-to-use bead-based platform for ultra-sensitive and low cost detection of analytes. Now, you can quantify multiple analytes in a single well using our new AlphaPlex™ technology - for faster and more accurate results, even when using low sample volumes. AlphaPlex is a homogeneous, all-in-one-well multiplexing technology that provides highly sensitive detection of a wide range of analytes in a simple no-wash assay format.

The EnVision® Multilabel Plate Reader delivers the ultimate in performance and reliability across all detection technologies, including Alpha Technology. If you already have an EnVision system then find out how easy it can be to enable the next step in Alpha Technology.

	EnVision Xcite or EnVision HTS with Standard Alpha <i>Product #2102* onwards</i>		EnVision Xcite or EnVision HTS without Standard Alpha <i>Product #2102* onwards</i>	
	Single Detector**	Dual Detector (HTS)***	Single Detector**	Dual Detector (HTS)***
	REQUIRED ACCESSORIES			
EnVision Alpha Standard Field Kit #2102-8330	Not needed	Not needed	Needed	Needed
AlphaPlex Filters Resorufine/Amplex Red FP 535/40 #2100-5570 Europium emission #2100-5090	Needed	Needed	Needed	Needed
Dual Alpha Mirror Module #2102-5900	Not needed	Needed for simultaneous measurement	Not needed	Needed for simultaneous measurement
EnVision Software Upgrade Kit**** #2101-8020 or EnVision Computer Upgrade Kit (Including Latest Software) #2104-8040	Software version 1.13 needs to be installed. Where older versions are currently installed, an upgrade will be needed.	Software version 1.13 needs to be installed. Where older versions are currently installed, an upgrade will be needed.	Software version 1.13 needs to be installed. Where older versions are currently installed, an upgrade will be needed.	Software version 1.13 needs to be installed. Where older versions are currently installed, an upgrade will be needed.

* We no longer offer service for product numbers #2102 and lower. Please be advised that although you can still upgrade, you do so at your own risk.

** Sequential read only

*** Sequential and dual read

**** If your computer was purchased before 2012, please check with your PerkinElmer representative as to whether you will require a new computer.

AlphaPlex is currently not available for the EnVision Xcite or EnVision HTS with Turbo Alpha, or for the EnSpire® Multimode Plate Reader or EnSight™ Multimode Plate Reader.

PerkinElmer, Inc.
940 Winter Street
Waltham, MA 02451 USA
P: (800) 762-4000 or
(+1) 203-925-4602
www.perkinelmer.com



For a complete listing of our global offices, visit www.perkinelmer.com/ContactUs

Copyright © 2014 PerkinElmer, Inc. All rights reserved. PerkinElmer® is a registered trademark of PerkinElmer, Inc. All other trademarks are the property of their respective owners.